


I. Barocycler Setup and Operation

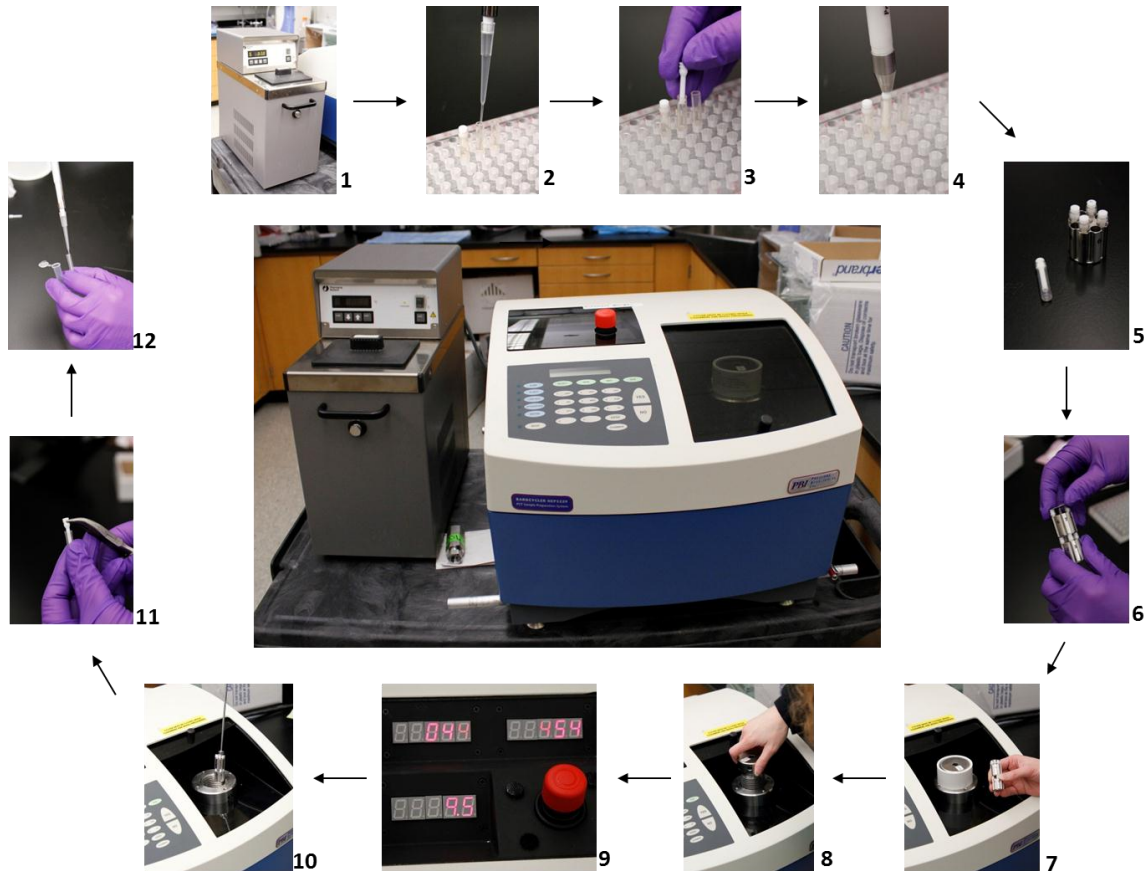

Barocycler Set-up and Operation (abridged from PBI User Manual for NEP 3229)

1. Turning on chiller to equilibrate at 37°C
2. Loading the PCT micro-tubes with gel loading pipette tips
3. Placing the cap into the PCT microtube
4. Capping with the high-pressure cap-insert tool
5. Loading and balancing the carousel
6. Assembling the carousels
7. Loading the carousel assembly
8. Capping the high pressure chamber
9. Running a cycle
10. Removing the carousel assembly with the magnetic wand
11. Removing the cap with vendor supplied pliers
12. Transferring sample with a gel loading pipette tip